



Contec ProChlor SATWIPE System

Meltblown polypropylene wipes with stabilised hypochlorous acid blended with purified water

Contec ProChlor is a revolutionary new sporicide achieving a 100% kill against spores in under 1 min. It is validated for sporicidal transfer disinfection in practice-mimicking conditions, achieving a log 5 kill on syringe paper samples and complete kill on plastic syringe samples, in 1 min.

A blend of hypochlorous acid in purified water, Contec ProChlor is efficacious against bacteria, fungi, moulds, yeasts, viruses and spores.

Contec ProChlor is provided as a point-of-use SATWIPE system. Each double bag pack contains a canister of 120 x 34gsm nonsterile meltblown polypropylene wipes and 700ml of either sterile or filtered Contec ProChlor, in a 1L bottle.

Simply add 700ml of ProChlor to the dry roll, 5 mins before use, and use the presaturated wipes within 8 hours.

Contec Sterile ProChlor is aseptic filled under UDAF grade A in a grade B cleanroom into pre-irradiated components. Contec ProChlor is 0.2 micron filtered and filled in a Grade B/C environment. The canister of dry wipes is packed in a Grade C cleanroom.



FEATURES	BENEFITS
Impregnated wipes at point-of-use	• Meets requirements of Aseptic Transfer Guidelines
Log 5 kill against <i>Bacillus subtilis</i> in 60 seconds	• Fast acting so suitable for transfer disinfection
120 polypropylene wipes in a canister	• Reduces amount of packaging generated • Fast and easy to use • High quality wipe for transfer disinfection
Contains no quaternary ammonium or surfactant	• Very low residue, saving time on residue removal
No hazard classification	• Good operator acceptability
Double bagged in linear tear packaging	• Each bag is easy to open even when wearing gloves



Product Specification

Contec SATWIPES with Sterile ProChlor

700ml capped - 2000ppm hypochlorous acid in purified water

Ready to use with a 2 year shelf life

Aseptic filled under UDAF grade A in a grade B cleanroom

34gsm polyester / cellulose wipes 150 x 200mm

Packed in a grade C cleanroom

Ready to use with a 2 year shelf life

Double bagged in polyethylene linear tear bags

Relevant certificates of analysis and sterility are provided with every batch. Full supporting documentation information is available on request.

Contec SATWIPES with ProChlor

700ml capped - 2000ppm hypochlorous acid in purified water

Ready to use with a 2 year shelf life

0.2 micron filtered in a grade B / C cleanroom

34gsm polyester / cellulose wipes 150 x 200mm

Packed in a grade C cleanroom

Ready to use with a 2 year shelf life

Double bagged in polyethylene linear tear bags

Efficacy

Contec ProChlor has been shown to have good efficacy against spores using standard laboratory tests. Contec ProChlor SATWIPES have also been tested against the EN 16615 test for presaturated wipes against bacteria, fungi and bacterial spores.

Test	Description	Log Reduction	Time	Test	Description	Log Reduction	Time
EN1276	<i>E. hirae</i>	> log 7	1 min	EN13704	<i>B. subtilis</i>	> log 6	1 min
EN1276	<i>S. aureus</i>	> log 7	1 min	EN13704	<i>B. pumilis</i>	>3.88*	1 min
EN1276	<i>P. aeruginosa</i>	> log 7	1 min	EN13704	<i>B. cereus</i>	>3.24*	1 min
EN1276	<i>E. coli</i>	>log 7	1 min	EN13697	<i>B. subtilis</i>	> log 6	1 min
EN14476	Vaccinia virus	log 5	30 secs	EN16615	<i>S. aureus</i>	7.24 test area 1**	1 min
EN14476	Adenovirus	4.67	30 secs	EN16615	<i>P. aeruginosa</i>	7.45 test area 1**	1 min
EN1650	<i>A. niger</i>	> log 4	1 min	EN16615	<i>A. brasiliensis</i>	6.13 test area 1**	1 min
EN1650	<i>C. albicans</i>	> log 4	1 min	EN16615	<i>A. subtilis</i>	7.11 test area 1**	1 min

* Complete kill ** Test fields 2-4 average of 0 cfu.

Product Application

Pour 700ml Contec ProChlor into the centre of the wipe roll. Close lid and invert canister multiple times for a 20 second period. Wait 5 minutes this will ensure saturation of all wipes. Following the 5-minute saturation time, some liquid will remain in the canister, this is intentional to ensure efficient wetting of the wipes. Once saturated with ProChlor, the wipes are suitable for sporocidal transfer disinfection. Wipes should be used within 8 hours & immediately disposed of. Dispose of remaining fluid and canister according to facility procedures.

Part No.	Description	Size	Packaging
XSWPP02PC	Contec Meltblown Polypropylene SATWIPES System For use with Contec Sterile ProChlor	150 x 200mm	120 wipes per canister; 10 canisters per case
FSWPP02PC	Contec Meltblown Polypropylene SATWIPES System For use with Contec ProChlor	150 x 200mm	120 wipes per canister; 10 canisters per case

Use biocides safely. Always read the label and product information before use.